Package 'topdownr'

April 12, 2018

Title Investigation of Fragmentation Conditions in Top-Down Proteomics

Version 1.0.0

Description The topdownr package allows automatic and systemic investigation of fragment conditions. It creates Thermo Orbitrap Fusion Lumos method files to test hundreds of fragmentation conditions. Additionally it provides functions to analyse and process the generated MS data and determine the best conditions to maximise overall fragment coverage.

Depends R (>= 3.4), methods, BiocGenerics (>= 0.20.0), Biostrings (>= 2.42.1), S4Vectors (>= 0.12.2)

- **Imports** stats, tools, utils, Biobase, Matrix (>= 1.2.10), MSnbase (>= 2.3.10), ggplot2 (>= 2.2.1), mzR (>= 2.11.4)
- Suggests topdownrdata (>= 0.2), knitr, ranger, testthat, BiocStyle

License GPL (>= 3)

URL https://github.com/sgibb/topdownr/

BugReports https://github.com/sgibb/topdownr/issues/

LazyData true

VignetteBuilder knitr

Roxygen list(markdown=TRUE)

RoxygenNote 6.0.1.9000

biocViews Infrastructure, Proteomics, MassSpectrometry, Coverage

Encoding UTF-8

NeedsCompilation no

Author Sebastian Gibb [aut, cre] (0000-0001-7406-4443), Pavel Shliaha [aut], Ole Nørregaard Jensen [aut]

Maintainer Sebastian Gibb <mail@sebastiangibb.de>

R topics documented:

topdownr-package					•	•				•								•					2
AbstractTopDownSet-class										•							•						3
createTngFusionMethFiles							•			•									 				6
defaultMs1Settings	•	•		•			•	•		•	•	•		•	•	•	•		•				8

topdownr-package

FragmentViews-class	. 8
NCBSet-class	. 10
readTopDownFiles	. 12
tds	. 14
TopDownSet-class	. 14
writeMethodXmls	. 18
	20

Index

topdownr-package Investigation of Fragmentation Conditions in Top-Down Proteomics

Description

The topdownr package allows automatic and systemic investigation of fragment conditions. It creates Thermo Orbitrap Fusion Lumos method files to test hundreds of fragmentation conditions. Additionally it provides functions to analyse and process the generated MS data and determine the best conditions to maximise overall fragment coverage.

Details

The usage of the topdownr package is demonstrated in the following vignettes:

- Generate .meth files prior data acquisition for the Thermo Orbitrap Fusion Lumos MS devise: vignette("data-generation", package="topdownr").
- How to analyse top-down fragmenation data: vignette("analysis", package="topdownr")

Author(s)

Sebastian Gibb <mail@sebastiangibb.de>, Pavel Shliaha <pavels@bmb.sdu.dk>, Ole Nørregaard Jensen <jenseno@bmb.sdu.dk>

References

https://github.com/sgibb/topdownr/

See Also

Useful links:

- https://github.com/sgibb/topdownr/
- Report bugs at https://github.com/sgibb/topdownr/issues/

AbstractTopDownSet-class

rowViews(object, ...)

The AbstractTopDownSet class

Description

Abstract/VIRTUAL parent class for TopDownSet and NCBSet to provide common interface.

Usage

```
## S4 method for signature 'AbstractTopDownSet,ANY,ANY,ANY'
x[i, j, ..., drop = FALSE]
## S4 method for signature 'AbstractTopDownSet,ANY,missing'
x[[i, j, ...]]
## S4 replacement method for signature 'AbstractTopDownSet,ANY,missing'
x[[i, j, ...]] <- value
## S4 method for signature 'AbstractTopDownSet'
x$name
## S4 replacement method for signature 'AbstractTopDownSet'
x$name <- value
## S4 method for signature 'AbstractTopDownSet'
assayData(object)
## S4 method for signature 'AbstractTopDownSet'
colData(object)
## S4 replacement method for signature 'AbstractTopDownSet'
colData(object, ...) <- value</pre>
## S4 method for signature 'AbstractTopDownSet'
conditionData(object, ...)
## S4 replacement method for signature 'AbstractTopDownSet'
conditionData(object, ...) <- value</pre>
## S4 method for signature 'AbstractTopDownSet'
dim(x)
## S4 method for signature 'AbstractTopDownSet'
dimnames(x)
## S4 method for signature 'AbstractTopDownSet'
removeEmptyConditions(object)
## S4 method for signature 'AbstractTopDownSet'
```

```
## S4 method for signature 'AbstractTopDownSet'
show(object)
## S4 method for signature 'AbstractTopDownSet'
summary(object, what = c("rows", "columns"),
    ...)
```

Arguments

i, j	numeric, logical or character, indices specifying elements to extract or replace.
drop	logical, currently ignored.
value	replacment value.
name	character name of an (non)existing column in colData.
object, x	AbstractTopDownSet
what	character, specifies whether "rows" or "columns" should be summarized.
	arguments passed to internal/other methods.

Details

This class just provides a common interface. It is not intended for direct use by the user. Please see TopDownSet for an example usage of its child class.

Value

This is an *Abstract/VIRTUAL* class to provide a common interface for TopDownSet and NCBSet. It is not possible to create an AbstractTopDownSet object.

Methods (by generic)

• [: Subset operator.

For i numeric, logical or character vectors or empty (missing) or NULL are supported. Subsetting is done on the fragment/bond (row) level. character indices could be names (e.g. c("a1", "b1", "c1", "c2", "c3")) or types (e.g. c("c", "x")) of the fragments for TopDownSet objects, or names of the bonds (e.g. c("bond001")) for NCBSet objects. j could be a numeric or logical vector and subsetting is done on the condition/run (column) level.

• [[: Subset operator.

i could be a numeric or logical vector and subsetting is done on the condition/run (column) level.

- [[<-: Setter for a column in the colData slot.
 The [[< operator is used to add/replace a single column of the colData DataFrame.
- \$: Accessor for columns in the colData slot.

The \$ simplifies the accession of a single column of the colData. It is identical to the [[operator.

• \$<-: Setter for a column in the colData slot.

The \$< - operator is used to add/replace a single column of the colData DataFrame. It is identical to the [[< - operator.

```
4
```

assayData: Accessor for the assay slot.

Returns a Matrix::dgCMatrix that stores the intensity/coverage information of AbstractTop-DownSet object.

• colData: Accessor for the colData slot.

Returns a S4Vectors::DataFrame that stores metadata for the conditons/runs (columns) of the AbstractTopDownSet object.

- colData<-: Setter for the colData slot.
 Replaces metadata for the conditons/runs (columns) of the AbstractTopDownSet object.
- conditionData: Accessor for the colData slot. An alias for colData.
- conditionData<-: Setter for the colData slot. An alias for colData<-.
- dim: Accessor for dimensions. Returns a numeric with number of fragments/bonds (rows) and conditions/runs (columns).
- dimnames: Accessor for dimension names.
 Returns a list with names for the fragments/bonds (rows) and for the conditions/runs (columns).
- removeEmptyConditions: Remove empty conditions/runs.
 Removes conditions/runs (columns) without any intensity/coverage information from the AbstractTopDownSet object. It returns a modified AbstractTopDownSet object.
- rowViews: Accessor for the rowViews slot.
 Depending on the implementation it returns an FragmentViews object for TopDownSet objects or an Biostrings::XStringViews object for NCBSet objects.
- summary: Summary statistics.

Returns a matrix with some statistics: number of fragments, total/min/first quartile/median/mean/third quartile/maximum of intensity values.

Slots

- rowViews Biostrings::XStringViews, information about fragments/bonds (name, type, sequence, mass, charge), see Biostrings::XStringViews and FragmentViews for details.
- colData S4Vectors::DataFrame, information about the MS2 experiments and the fragmentation conditions.
- assay Matrix::dgCMatrix, intensity/coverage values of the fragments/bonds. The rows correspond to the fragments/bonds and the columns to the condition/run. It just stores values that are different from zero.
- files character, files that were imported.
- tolerance double, tolerance in *ppm* that were used for matching the experimental m/z values to the theoretical fragments.

processing character, log messages.

Author(s)

Sebastian Gibb <mail@sebastiangibb.de>

See Also

- TopDownSet and NCBSet which both implement/use this interface. These manual pages also provide some example code.
- FragmentViews (and Biostrings::XStringViews) for the row view interface.
- Matrix::dgCMatrix for technical details about the intensity/coverage storage.

Examples

```
## Because AbstractTopDownSet is a VIRTUAL class we could not create any
## object of it. Here we demonstrate the usage with an TopDownSet that
## implements the AbstractTopDownSet interface. See `?"TopDownSet-class"` for
## more details an further examples.
## Example data
data(tds, package="topdownr")
tds
head(summary(tds))
# Accessing slots
rowViews(tds)
colData(tds)
assayData(tds)
# Accessing colData
tds$Mz
tds$FilterString
# Subsetting
# First 100 fragments
tds[1:100]
# All c fragments
tds["c"]
# Just c 152
tds["c152"]
# Condition 1 to 10
tds[, 1:10]
```

createTngFusionMethFiles

Windows specific functions.

Description

The functions runXmlMethodChanger and runScanHeadsman call XmlMethodChanger.exe and ScanHeadsman.exe with the corresponding arguments. The only work on Windows (maybe on Linux + wine as well but that was never tested).

Usage

```
createTngFusionMethFiles(template, xml = list.files(pattern = ".*\\.xml$"),
    executable = "XmlMethodChanger.exe", verbose = interactive())
runXmlMethodChanger(template, xml = list.files(pattern = ".*\\.xml$"),
```

```
executable = "XmlMethodChanger.exe", verbose = interactive())
```

```
runScanHeadsman(path = ".", executable = "ScanHeadsman.exe")
```

Arguments

template	character, path to template .meth file.
xml	character, vector of path to .xml files.
executable	character, path to the XmlMethodChanger.exe or ScanHeadsman.exe exe- cutable.
verbose	logical, if TRUE a progress bar is shown.
path	character, path to the directory containing the .raw files.

Details

runXmlMethodChanger applies 'XmlMethodChanger.exe' on all given XML files generated with writeMethodXmls() to create .meth files from a template.

runScanHeadsman calls ScanHeadsman.exe on a given directory containing .raw files. ScanHeadsman.exe extracts the method and scan header data into .experiments.csv and .txt files, respectively.

Value

Nothing. Used for its side effects.

References

XmlMethodChanger source code: https://github.com/thermofisherlsms/meth-modifications/ ScanHeadsman source code: https://bitbucket.org/caetera/scanheadsman

See Also

writeMethodXmls()

Examples

End(Not run)

defaultMs1Settings Settings for MS1/2 parameters.

Description

This functions create the default settings for writeMethodXmls().

Usage

```
defaultMs1Settings(...)
```

```
defaultMs2Settings(...)
```

Arguments

...

named arguments that should be overwritten.

Value

A named list of settings.

Examples

```
# all default MS1 settings
defaultMs1Settings()
# overwrite FirstMass and set AgcTarget
```

```
defaultMs1Settings(FirstMass=100, AgcTarget=c(1e4, 1e5))
```

```
# all default MS2 settings
defaultMs2Settings()
```

```
# overwrite AgcTarget
defaultMs2Settings(AgcTarget=c(1e4, 1e5))
```

FragmentViews-class The FragmentViews class

Description

The FragmentViews class is a basic container for storing a set of views (start/end locations) on the same peptides/protein sequence. Additionally it keeps information about mass, type and charge of the fragments.

Usage

```
FragmentViews(sequence, mass, type, z = 1L, start = NULL, end = NULL,
width = NULL, names = NULL, metadata = list())
## S4 method for signature 'FragmentViews'
show(object)
```

Arguments

sequence	character/ Biostrings:: AAString, complete protein/peptide sequence.
mass	double, mass of the fragments, same length as start/end/width.
type	character, type of the fragments, same length as start/end/width'.
Z	integer, charge of the fragments, length one or same length as start/end/width'.
start	integer, start positions of the fragments. At least two of start/end/width' has to be given.
end	integer, end positions of the fragments. At least two of start/end/width' has to be given.
width	integer, width positions of the fragments. At least two of start/end/width' has to be given.
names	character, names of the fragments, same length as start/end/width'.
metadata	list, metadata like modifications.
object	FragmentViews

Details

FragmentViews extends Biostrings::XStringViews. In short it combines an IRanges::IRanges object to store start/end location on a sequence, an Biostrings::AAString object.

Value

An FragmentViews object.

Functions

• FragmentViews: Constructor In general it is not necessary to call the constructor manually. See readTopDownFiles() instead.

Coercion

as(object, "data.frame"): Coerce an FragmentViews object into a data.frame.

Author(s)

Sebastian Gibb <mail@sebastiangibb.de>

See Also

Biostrings::XStringViews

Examples

NCBSet-class

Description

The NCBSet class is a container for a top-down proteomics experiment similar to the TopDownSet but instead of intensity values it just stores the information if a bond is covered by a N-terminal (encoded as 1), a C-terminal (encoded as 2) and/or bidirectional fragments (encoded as 3).

Usage

```
## S4 method for signature 'NCBSet'
bestConditions(object, n = ncol(object), minN = 0L, ...)
## S4 method for signature 'NCBSet'
fragmentationMap(object, nCombinations = 10,
    cumCoverage = TRUE, labels = colnames(object), ...)
## S4 method for signature 'NCBSet'
show(object)
## S4 method for signature 'NCBSet'
summary(object, what = c("conditions", "bonds"), ...)
```

Arguments

object	NCBSet
n	integer, max number of combinations/iterations.
minN	integer, stop if there are less than minN additional fragments
nCombinations	integer, number of combinations to show (0 to avoid plotting them at all).
cumCoverage	logical, if TRUE (default) cumulative coverage of combinations is shown.
labels	character, overwrite x-axis labels.
what	character, specifies whether "conditions" (columns; default) or "bonds" (rows) should be summarized.
	arguments passed to internal/other methods. added. newly added fragments third column: number of newly covered bonds.

Value

An NCBSet object.

Methods (by generic)

• bestConditions: Best combination of conditions.

Finds the best combination of conditions for highest coverage of bonds. Use n to limit the number of iterations and combinations that should be returned. If minN is set at least minN fragments have to be added to the combinations. The function returns a 2-column matrix. The first column contains the index of the condition (column number) and the second one the newly added number of fragments.

NCBSet-class

• fragmentationMap: Plot fragmentation map.

Plots a fragmentation map of the Protein. Use nCombinations to add another plot with nCombinations combined conditions. If cumCoverage is TRUE (default) these combinations increase the coverage cumulatively.

• summary: Summary statistics.

Returns a matrix with some statistics: number of fragments, total/min/first quartile/median/mean/third quartile/maximum of intensity values.

Slots

- rowViews Biostrings::XStringViews, information about bonds (name, start, end, width, sequence), see Biostrings::XStringViews for details.
- colData S4Vectors::DataFrame, information about the MS2 experiments and the fragmentation conditions.
- assay Matrix::dgCMatrix, coverage values of the bonds. The rows correspond to the bonds and the columns to the condition/run. It just stores values that are different from zero. If a bond is covered by an N-terminal fragment its encoded with 1, by an C-terminal fragmentl with 2 and by both fragment types/bidirectional by 3 respectively.
- files character, files that were imported.
- tolerance double, tolerance in *ppm* that were used for matching the experimental mz values to the theoretical fragments.

processing character, log messages.

Author(s)

Sebastian Gibb <mail@sebastiangibb.de>

See Also

- An NCBSet is generated from an TopDownSet object.
- Biostrings::XStringViews for the row view interface.
- Matrix::dgCMatrix for technical details about the coverage storage.

Examples

```
## Example data
data(tds, package="topdownr")
```

Aggregate technical replicates
tds <- aggregate(tds)</pre>

Coercion into an NCBSet object
ncb <- as(tds, "NCBSet")</pre>

ncb

head(summary(ncb))

Accessing slots
rowViews(ncb)
colData(ncb)
assayData(ncb)

```
# Accessing colData
ncb$Mz
# Subsetting
# First 100 bonds
ncb[1:100]
# Just bond 152
ncb["bond152"]
# Condition 1 to 10
ncb[, 1:10]
# Plot fragmentation map
fragmentationMap(ncb)
```

readTopDownFiles Read top-down files.

Description

It creates an TopDownSet object and is its only constructor.

Usage

```
readTopDownFiles(path, pattern = ".*", type = c("a", "b", "c", "x", "y",
    "z"), modifications = c("Carbamidomethyl", "Acetyl", "Met-loss"),
    adducts = data.frame(), neutralLoss = MSnbase::defaultNeutralLoss(),
    sequenceOrder = c("original", "random", "inverse"), tolerance = 5e-06,
    dropNonInformativeColumns = TRUE, sampleColumns = c("Mz", "AgcTarget",
    "EtdReagentTarget", "EtdActivation", "CidActivation", "HcdActivation"),
    verbose = interactive())
```

Arguments

path	character, path to directory that contains the top-down files.
pattern	character, a filename pattern, the default .* means all files.
type	character, type of fragments, currently $a - c$ and $x - z$ are supported, see MSnbase::calculateFragmer for details.
modifications	character, unimod names of modifications that should be applied. Currenlty just <i>Acetyl</i> (Unimod:1 but just protein N-term), <i>Carbamidomethyl</i> (Unimod:4) and <i>Met-loss</i> (Unimod:765) are supported. <i>Met-loss</i> removes M (if followed by A, C, G, P, S, T, or V; (see also http://www.unimod.org/modifications_view.php?editid1=1, http://www.unimod.org/modifications_view.php?editid1=4, and http://www.unimod.org/modification for details)). Use NULL to disable all modifications.
adducts	data.frame, with 3 columns, namely: mass, name, to, see details section.
neutralLoss	list, neutral loss that should be applied, see MSnbase::calculateFragments() and MSnbase::defaultNeutralLoss() for details.

12

sequenceOrder	character, order of the sequence before fragment calculation and matching is done. "original" doesn't change anything. "inverse" reverse the sequence and "random" arranges the amino acid sequence at random.				
tolerance	double, tolerance in <i>ppm</i> that is used to match the theoretical fragments with				
the observed ones. dropNonInformativeColumns					
	logical, should columns with just one identical value across all runs be removed?				
sampleColumns	character, column names of the colData() used to define a sample (technical replicate). This is used to add the Sample column (used for easier aggregation, etc.).				
verbose	logical, verbose output?				

Details

readTopDownFiles reads and processes all top-down files, namely:

- .fasta (protein sequence)
- .mzML (spectra)
- .experiments.csv (method/fragmentation conditions)
- .txt (scan header information)

adducts: *Thermo's Xtract* allows some mistakes in deisotoping, mostly it allows +/- C13-C12 and +/- H+. The adducts argument takes a data.frame with the mass to add, the name that should assign to these new fragments and an information to whom the modification should be applied, e.g. for H+ on z, data.frame(mass=1.008, name="zpH", to="z").

Please note: The adducts are added to the output of MSnbase::calculateFragments(). That has some limitations, e.g. neutral loss calculation could not be done in topdownr-package. If neutral loss should be applied on adducts you have to create additional rows, e.g.: data.frame(mass=c(1.008, 1.008), name=c("

Value

A TopDownSet object.

See Also

MSnbase::calculateFragments(), MSnbase::defaultNeutralLoss()

Examples

```
if (require("topdownrdata")) {
    # add H+ to z and no neutral loss of water
    tds <- readTopDownFiles(
        topdownrdata::topDownDataPath("myoglobin"),
        ## Use an artifical pattern to load just the fasta
        ## file and files from m/z == 1211, ETD reagent
        ## target 1e6 and first replicate to keep runtime
        ## of the example short
        pattern=".*fasta.gz$|1211_.*1e6_1",
        adducts=data.frame(mass=1.008, name="zpH", to="z"),
        neutralLoss=MSnbase::defaultNeutralLoss(
            disableWaterLoss=c("Cterm", "D", "E", "S", "T")),
        tolerance=25e-6
    )
}</pre>
```

tds

Description

An example data set for topdownr. It is just a subset of the myoglobin dataset available in topdownrdata::topdownrdata-package.

Usage

tds

Format

A TopDownSet with 2700 fragments and 351 conditions.

Details

It was created as follows:

```
tds <- readTopDownFiles(
   topdownrdata::topDownDataPath("myoglobin"),
   ## Use an artifical pattern to load just the fasta
   ## file and files from m/z == 1211, ETD reagent
   ## target 1e6 and first replicate to keep runtime
   ## of the example short
   pattern=".*fasta.gz$|1211_.*1e\+06_1",
   adducts=data.frame(mass=1.008, name="zpH", to="z"),
   neutralLoss=MSnbase::defaultNeutralLoss(
        disableWaterLoss=c("Cterm", "D", "E", "S", "T")),
   tolerance=25e-6)</pre>
```

Source

Subset taken from the topdownrdata::topdownrdata-package package.

TopDownSet-class The TopDownSet class

Description

The TopDownSet class is a container for a whole top-down proteomics experiment.

TopDownSet-class

Usage

```
## S4 method for signature 'TopDownSet'
aggregate(x, by = x$Sample, ...)
## S4 method for signature 'TopDownSet'
filterCv(object, threshold, by = object$Sample, ...)
## S4 method for signature 'TopDownSet'
filterInjectionTime(object, maxDeviation = log2(3),
 keepTopN = 2, by = object$Sample, ...)
## S4 method for signature 'TopDownSet'
filterIntensity(object, threshold, relative = TRUE,
  ...)
## S4 method for signature 'TopDownSet'
filterNonReplicatedFragments(object, minN = 2,
 by = object$Sample, ...)
## S4 method for signature 'TopDownSet'
normalize(object, method = "TIC", ...)
## S4 method for signature 'TopDownSet'
show(object)
## S4 method for signature 'TopDownSet'
summary(object, what = c("conditions", "fragments"),
  ...)
```

Arguments

x, object	TopDownSet
by	list, grouping variable, in general it refers to technical
threshold	double, threshold variable.
maxDeviation	double, maximal allowed deviation in the log2 injection time in ms in compar- ison to the median ion injection time.
keepTopN	integer, how many technical replicates should be kept?
relative	logical, if relative is TRUE all fragments with intensity below threshold * max(intensity) per fragment are removed, otherwise all fragments below threshold are removed.
minN	numeric, if less than minN of a fragment are found across technical replicates it is removed.
method	character, normalisation method, currently just "TIC" for Total Ion Current normalisation of the scans/conditions (column-wise normalisation) is supported.
what	character, specifies whether "conditions" (columns; default) or "fragments" (rows) should be summarized.
	arguments passed to internal/other methods. replicates (that's why the default is the "Sample" column in colData).

Details

See vignette("analysis", package="topdownr") for a detailed example how to work with TopDownSet objects.

Value

An TopDownSet object.

Methods (by generic)

• aggregate: Aggregate conditions/runs.

Aggregates conditions/runs (columns) in an TopDownSet object by a user-given value (default is the "Sample" column of colData which has the same value for technical replicates). It combines intensity values and numeric metadata of the grouped conditions/runs (columns) by mean and returns a reduced TopDownSet object.

• filterCv: Filter by CV.

Filtering is done by coefficient of variation across technical replicates (defined by the by argument). All fragments below a given threshold are removed. The threshold is the maximal allowed CV in percent (sd/mean * 100 < threshold).

• filterInjectionTime: Filter by ion injection time.

Filtering is done by maximal allowed deviation and just the technical keepTopN replicates with the lowest deviation from the median ion injection time are kept.

- filterIntensity: Filter by intensity.
 Filtering is done by removing all fragments that are below a given (absolute/relative) intensity threshold.
- filterNonReplicatedFragments: Filter by non-replicated fragments.

Filtering is done by removing all fragments that don't replicate across technical replicates.

• normalize: Normalise.

Applies *T*otal *I*on *C*urrent normalisation to a TopDownSet object. The normalisation ist done per scans/conditions (column-wise normalisation).

• summary: Summary statistics.

Returns a matrix with some statistics: number of fragments, total/min/first quartile/median/mean/third quartile/maximum of intensity values.

Slots

- rowViews FragmentViews, information about fragments (name, type, sequence, mass, charge), see FragmentViews for details.
- colData S4Vectors::DataFrame, information about the MS2 experiments and the fragmentation conditions.
- assay Matrix::dgCMatrix, intensity values of the fragments. The rows correspond to the fragments and the columns to the condition/run. It just stores values that are different from zero.
- files character, files that were imported.
- tolerance double, tolerance in *ppm* that were used for matching the experimental mz values to the theoretical fragments.

processing character, log messages.

TopDownSet-class

Coercion

'as(object, "MSnSet"): Coerce an TopDownSet object into an MSnbase::MSnSet object.'as(object, "NCBSet"): Coerce an TopDownSet object into an NCBSet object.

Author(s)

Sebastian Gibb <mail@sebastiangibb.de>

See Also

- FragmentViews for the row view interface.
- Matrix::dgCMatrix for technical details about the intensity storage.
- ?vignette("analysis", package="topdownr") for a full documented example of an analysis using topdownr.

Examples

```
## Example data
data(tds, package="topdownr")
```

tds

head(summary(tds))

Accessing slots
rowViews(tds)
colData(tds)
assayData(tds)

Accessing colData
tds\$Mz
tds\$FilterString

```
# Subsetting
```

First 100 fragments
tds[1:100]

All c fragments
tds["c"]

Just c 152 tds["c152"]

Condition 1 to 10 tds[, 1:10]

```
# Filtering
# Filter all intensities that don't have at least 10 % of the highest
# intensity per fragment.
tds <- filterIntensity(tds, threshold=0.1)</pre>
```

Filter all conditions with a CV above 30 % (across technical replicates)
tds <- filterCv(tds, threshold=30)</pre>

```
# Filter all conditions with a large deviation in injection time
tds <- filterInjectionTime(tds, maxDeviation=log2(3), keepTopN=2)
# Filter all conditions where fragments don't replicate
tds <- filterNonReplicatedFragments(tds)
# Normalise by TIC
tds <- normalize(tds)
# Aggregate technical replicates
tds <- aggregate(tds)
head(summary(tds))
# Coercion
as(tds, "NCBSet")
if (require("MSnbase")) {
    as(tds, "MSnSet")
}
```

writeMethodXmls Create Orbitrap Fusion method.xml files.

Description

This function is used to create Orbitrap Fusion method files from all combinations of a user-given set of MS1 and MS2 settings.

Usage

```
writeMethodXmls(ms1Settings, ms2Settings, groupBy = c("replication",
   "ETDReactionTime"), replications = 2, mz, massLabeling = TRUE,
   nMs2perMs1 = 10, duration = 0.5, randomise = TRUE,
   pattern = "method_%s.xml", verbose = interactive())
```

Arguments

ms1Settings	list, named list of MS1 settings, see e.g. defaultMs1Settings().
ms2Settings	list, named list of MS2 settings, see e.g. defaultMs2Settings().
groupBy	character, split files by groupBy entries.
replications	integer, number of replications of experiments.
mz	<pre>matrix, two columns (column 1: mass, column 2: z).</pre>
massLabeling	logical, should mz values used for ID labeling?
nMs2perMs1	integer, how many MS2 scans should be run after a MS1 scan?
duration	double, how long should the scan be?
randomise	logical, should the MS2 scan settings randomised?
pattern	character, file name pattern for the method.xml files.
verbose	logical, verbose output?

```
18
```

writeMethodXmls

Details

- ms1Settings: A list of MS1 settings. This has to be a named list. Valid MS1 settings are: c("FirstMass", "LastMass", "Microscans", "MaxITTimeInMS", "AgcTarget")
- ms2Settings: A list of MS2 settings. This has to be a named list. Valid MS2 settings are: c("ActivationType", "IsolationWindow", "EnableMultiplexIons", "EnableMSXIds", "MaxNoOfMultip
- groupBy: The groupBy parameter is used to split methods into different files. Valid entries are all settings that could be used in ms2Settings and "replication".
- massLabeling: The Orbitrap Fusion devices seems not to respect the start and end times of the runs given in the method.xml files. That's why it is nearly impossible to identify the run with its conditions based on the timings. If massLabeling is TRUE (default) the mass values given in mz are rounded to the first decimal place and the second to fourth decimal place is used as numeric identifier.
- pattern: The file name pattern used to name different method files. It must contain a "%s" that is replaced by the conditions defined in groupBy.

Value

An invisivble list with the MS1, MS2, runtimes and mz.

Author(s)

Sebastian Gibb <mail@sebastiangibb.de>, Pavel V. Shliaha <pavels@bmb.sdu.dk>

See Also

defaultMs1Settings(), defaultMs2Settings()

Examples

```
writeMethodXmls(defaultMs1Settings(FirstMass=400),
                defaultMs2Settings(),
                mz=cbind(mass=c(609.21, 700.45, 823.95), z=10),
                groupBy="ETDReactionTime",
                replications=1,
                pattern="method_firstmass_400_%s.xml")
```

Index

*Topic **datasets** tds, 14 *Topic package topdownr-package, 2[,AbstractTopDownSet,ANY,ANY,ANY-method (AbstractTopDownSet-class), 3 [[,AbstractTopDownSet,ANY,missing,-method (AbstractTopDownSet-class), 3 [[,AbstractTopDownSet,ANY,missing-method (AbstractTopDownSet-class), 3 [[<-,AbstractTopDownSet,ANY,missing,-method</pre> (AbstractTopDownSet-class), 3 [[<-,AbstractTopDownSet,ANY,missing-method</pre> (AbstractTopDownSet-class), 3 \$,AbstractTopDownSet-method (AbstractTopDownSet-class), 3 \$<-,AbstractTopDownSet-method</pre> (AbstractTopDownSet-class), 3

AbstractTopDownSet, 5 AbstractTopDownSet-class, 3 aggregate, TopDownSet-method (TopDownSet-class), 14 assayData, AbstractTopDownSet-method (AbstractTopDownSet-class), 3

coerce,FragmentViews,data.frame-method (FragmentViews-class), 8 coerce,TopDownSet,MSnSet-method (TopDownSet-class), 14 coerce,TopDownSet,NCBSet-method (TopDownSet-class), 14 colData(AbstractTopDownSet-class), 3 colData(), 13 colData,AbstractTopDownSet-method (AbstractTopDownSet-class), 3 colData<-(AbstractTopDownSet-class), 3</pre> defaultMs1Settings, 8 defaultMs1Settings(), 18, 19 defaultMs2Settings (defaultMs1Settings), 8 defaultMs2Settings(), 18, 19 dim,AbstractTopDownSet-method (AbstractTopDownSet-class), 3 dimnames, AbstractTopDownSet-method (AbstractTopDownSet-class), 3 filterCv (TopDownSet-class), 14 filterCv,TopDownSet-method (TopDownSet-class), 14 filterInjectionTime (TopDownSet-class), 14 filterInjectionTime,TopDownSet-method (TopDownSet-class), 14 filterIntensity (TopDownSet-class), 14 filterIntensity, TopDownSet-method (TopDownSet-class), 14 filterNonReplicatedFragments (TopDownSet-class), 14 filterNonReplicatedFragments,TopDownSet-method (TopDownSet-class), 14 fragmentationMap (NCBSet-class), 10 fragmentationMap,NCBSet-method (NCBSet-class), 10 FragmentViews, 5, 6, 9, 16, 17 FragmentViews (FragmentViews-class), 8 FragmentViews-class, 8

IRanges::IRanges, 9

INDEX

```
Matrix::dgCMatrix, 5, 6, 11, 16, 17
MSnbase::calculateFragments(), 12, 13
MSnbase::defaultNeutralLoss(), 12, 13
MSnbase::MSnSet, 17
```

NCBSet, 3-6, 10, 17 NCBSet-class, 10 normalize, TopDownSet-method (TopDownSet-class), 14

```
readTopDownFiles, 12
readTopDownFiles(), 9
removeEmptyConditions
    (AbstractTopDownSet-class), 3
removeEmptyConditions,AbstractTopDownSet-method
    (AbstractTopDownSet-class), 3
rowViews(AbstractTopDownSet-class), 3
rowViews,AbstractTopDownSet-class), 3
rowViews,AbstractTopDownSet-class), 3
runScanHeadsman
    (createTngFusionMethFiles), 6
runXmlMethodChanger
    (createTngFusionMethFiles), 6
```

```
tds, 14
topdownr (topdownr-package), 2
topdownr-package, 2, 13
topdownrdata::topdownrdata-package, 14
TopDownSet, 3–6, 10–12, 14, 16, 17
TopDownSet-class, 14
```

writeMethodXmls, 18
writeMethodXmls(), 7, 8